



LABCONCO CORPORATION 8811 Prospect Ave, Kansas City, MO 64132
(816) 333-8811 or (800) 821-5525 • (816) 363-0130 fax • labconco.com

Ceiling Enclosure Kit Installation

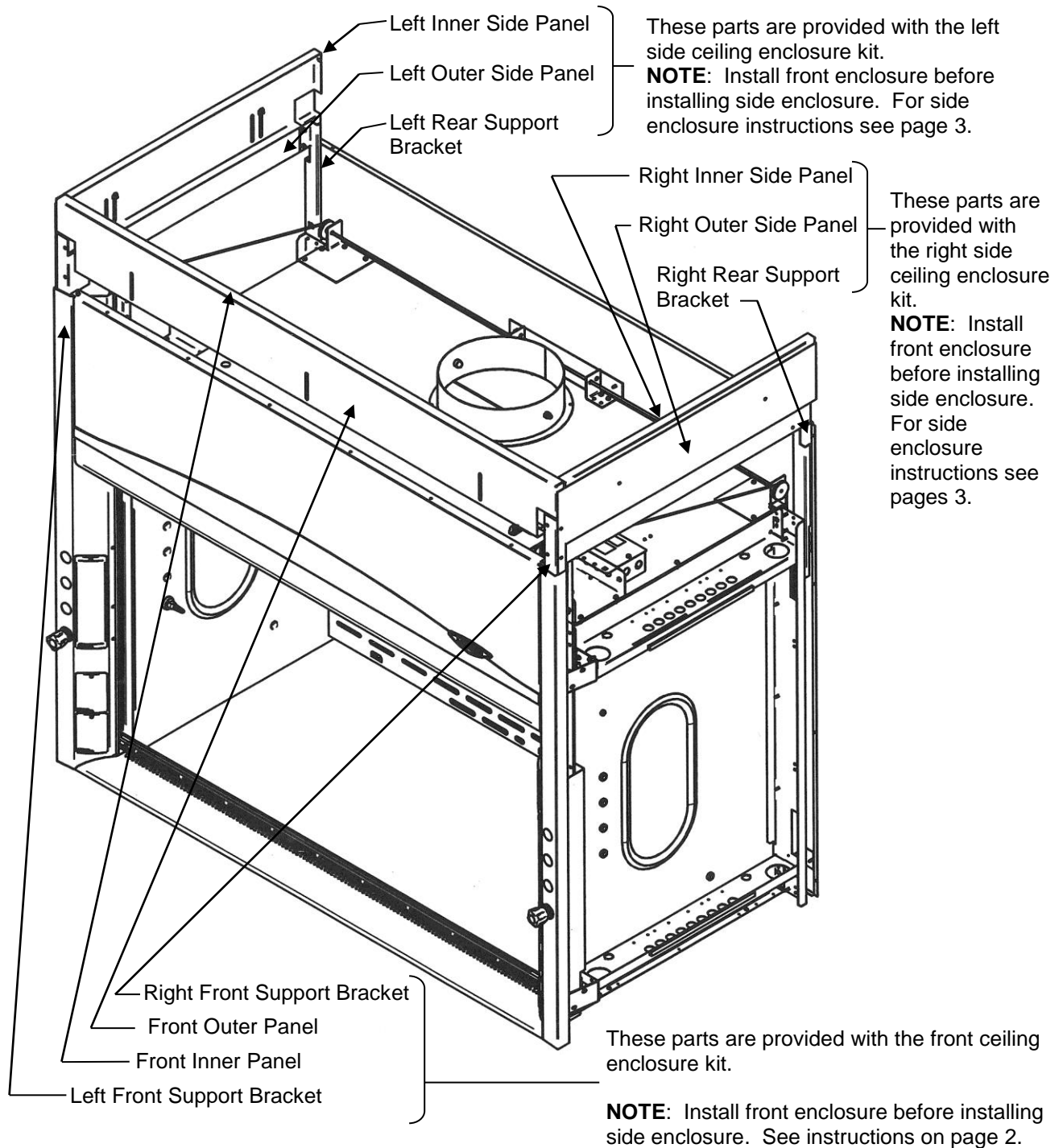

Front Panel Installation

1. Remove the front panel from the fume hood and lower sash.
2. Assemble the front inner panel to the front outer panel with flat washers and keps nuts, as shown.
3. Install left and right front support brackets to the ceiling enclosure front assembly using #10-24 keps nuts.
4. Drop assembly down into corner posts and fasten with #10-24 x .50 screws.

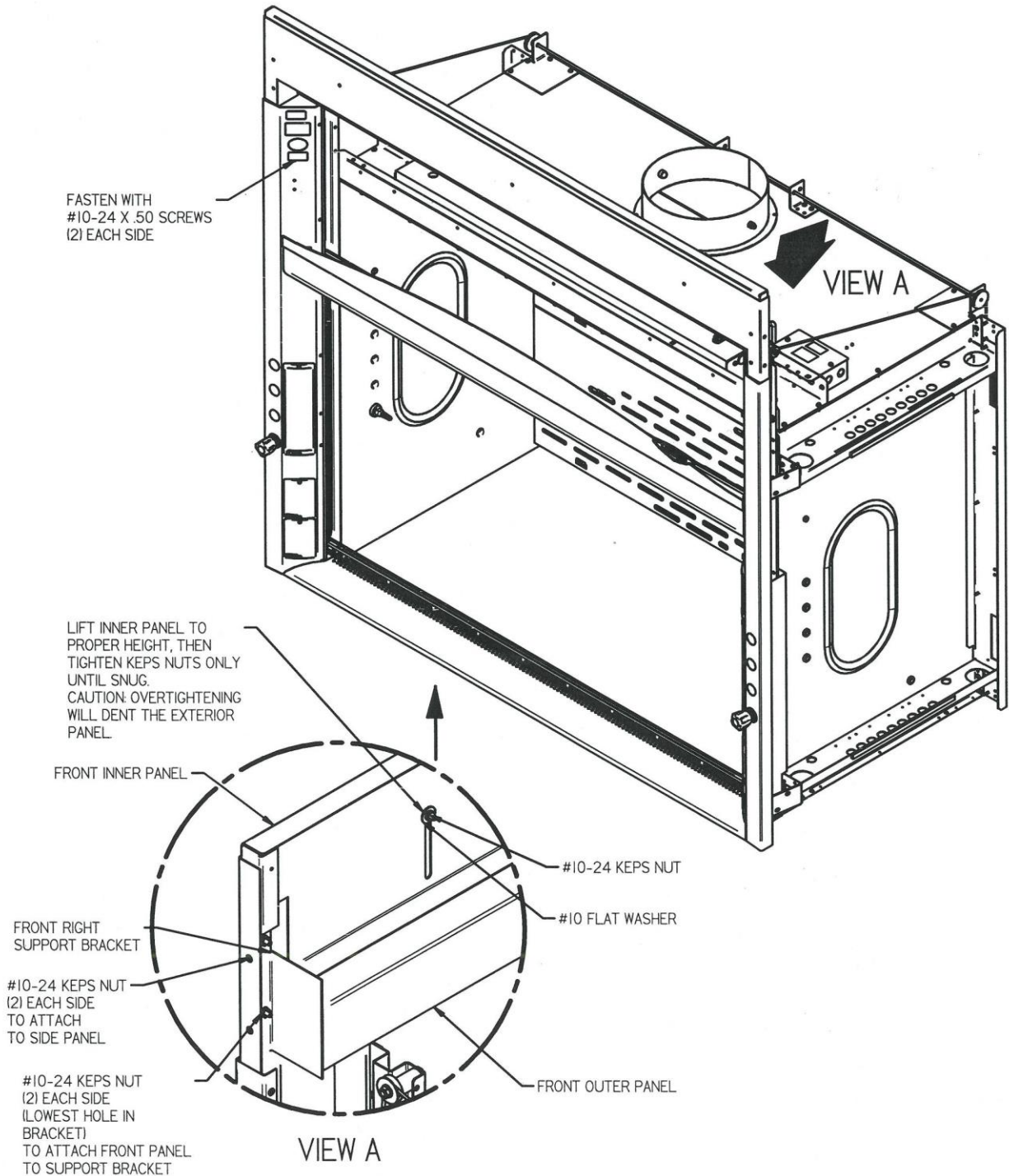

Side Panel Installation

NOTE: Install the front ceiling enclosure kit before installing the side ceiling enclosure kits.

The following instructions show the right side of fume hood. The left side enclosure kit installs in the same manner.

1. Remove the hood side panel.
2. Assemble the right rear support bracket to the right outer panel with (2) #10-24 keps nuts.
3. Assemble the inner side panel to the outer side panel with flat washers and keps nuts as shown below.
4. Assemble the side panel assembly (with inner panel pre-assembled to outer panel) and support bracket to the rear vertical channel and front panel assembly, using hardware as shown below.
5. Adjust the inner panel height to same height as front panel and tighten the two keps nuts. Secure with #10 x .50 Type AB screw.

